Informationsbroschüre des Herstellers für: Information Borchure supplied by Manufacturer:

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*1883: ECS GmbH Huettfeldstrasse 50, D-73430 Aalen

Contents

1. General Notes	2
2. Use, storage, maintenance	2
3. Cleaning and disinfection.....	2
4. Durability and spare parts	2
5. Area of application and protection, services and markings.....	2

1. General notes

The use of the face shield presupposes full knowledge of and compliance with this information brochure. The eye protection device is intended exclusively for the purposes stated in the information brochure.

EKASTU Safety GmbH assumes liability within the scope of its general terms and conditions. If you do not have them, they will be sent to you on request. The manufacturer is not liable for damage caused by failure to comply with the information brochure.

2. Use, Storage, Maintenance

The k1 face shield can be used for protection against mechanical impacts and for protection against liquids (drops and splashes e.g. acid and alkaline solutions). It can be worn in conjunction with half masks or filtering face pieces.

The visor of the k1 can be continuously tilted up. Make sure that the eye protection fits comfortably when you wear it. The visor is mounted on the holding device as follows: Insert the 3 screws that are included when ordering order No. 170 574, each through one of the pre-drilled holes in the holding device and through the corresponding recess on the top of the visor. Tighten all screws with supplied nut.

The visor should be stored and transported protected from direct sunlight, chemicals and scratching media.

3. Cleaning and disinfection

Clean the face shield with lukewarm water and addition of EKASTU cleaning fluid (0.5% solution) (order no. 266 962). After cleaning, rinse with clear water and dry with a soft cloth.

Disinfect the face shield with lukewarm water and addition of EKASTU disinfectant (2.0% solution) (order no. 266 942). After disinfection, rinse with clear water and dry with a soft cloth.

4. Durability and spare parts

Under normal conditions the visor of the face shield provides 1-2 years of adequate eye protection. For the model k1, replacement visors (order no. 170 604) can be ordered subsequently, which are replaced as described under point 2. Furthermore, screw replacement sets (order no. 140 090) can be ordered.

5. Area of application and protection, services and markings

The face shield is for your personal safety. EKASTU Safety GmbH guarantees the specified performance according to class and type. Please note that laboratory test values may differ considerably from those achieved in practice. The user must read and understand all instructions for use. Use the face shield exclusively in accordance with the information brochure, the applicable legal regulations and the safety regulations of the professional associations, in particular the rules of use according to DGUV rule 112-192 "Use of ocular protection and face protection".

The marking of the face protection device merely confirms that it complies with the basic requirements of the European standards EN 166 and is approved. That it is also approved for other purposes is indicated by letters and numbers on the lens as follows:

Visor's marking

Manufacturer	EKASTU Safety
Order number; Article description	170 604; k1
Certification mark	CE *
Number of the European standard	EN 166
Optical class	1
Identification symbol for field of use	3 (Liquids (droplets and splashes))
Identification symbol for mechanical strength / protection against mechanical impacts	S (increased robustness)

Holding devices's marking

Manufacturer	EKASTU Safety
Order number; Article description	170 574; k1
Certification mark	CE *
Number of the European standard	EN 166
Optical class	1
Identification symbol for field of use	3 (Liquids (droplets and splashes))
Identification symbol for mechanical strength / protection against mechanical impacts	S (increased robustness)

Attention:

The above-mentioned visor is shatterproof and scratch-resistant, but not unbreakable.

Check and clean it regularly. Scratched or cracked windows impair visibility and only provide a greatly reduced protective effect. They must therefore be replaced immediately. If you have any questions or doubts, contact your safety officer or the manufacturer.

In the case of persons with the appropriate disposition, allergic reactions can be triggered by contact of the skin with the wearer's body; in such cases a doctor should be consulted.

If the visor and holding device bear different markings for mechanical strength and protection against particles, the lower protection applies to the complete eye protection device.

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